



Maintenance for XLT U100

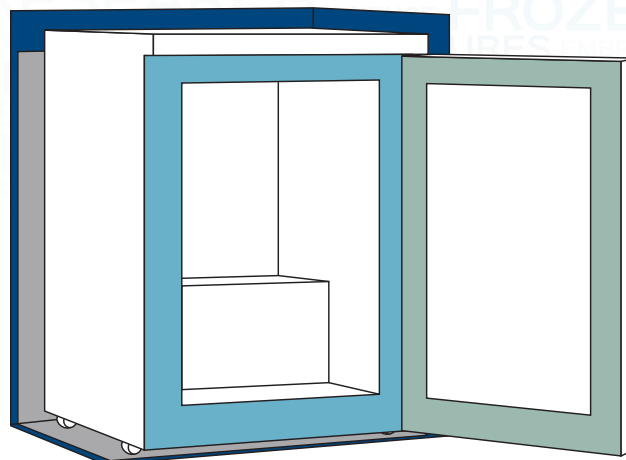

Maintenance										
Dark blue	Weekly	<p>Check placement:</p> <p>Min. 6 cm free space to the side of the freezer</p> <p>Min. 10 cm free space to the back of the freezer.</p> <p>If placed under a table, ensure that there is at least 6 cm free space above or a ventilation hole in the table or a gap between the table and the wall.</p>								
Light blue	Weekly	<p>Defrosting:</p> <p>Small amounts of loose ice can be removed using a plastic scraper.</p> <p>Defrosting has to be done regularly to prevent ice building</p> <p>The ice build depends on the use of the freezer.</p> <p>The more daily door openings - the more manual defrosting.</p> <table><tr><td>Door open/close per 24h</td><td>Advised monthly manual defrosting</td></tr><tr><td>< 5</td><td>1-2</td></tr><tr><td>< 10</td><td>2-4</td></tr><tr><td>< 15</td><td>4-6</td></tr></table>	Door open/close per 24h	Advised monthly manual defrosting	< 5	1-2	< 10	2-4	< 15	4-6
Door open/close per 24h	Advised monthly manual defrosting									
< 5	1-2									
< 10	2-4									
< 15	4-6									
Gray	Monthly	<p>Clean:</p> <p>The ventilation grille, the condenser, the compressor, the compressor compartment and the floor must also be kept free of dust and dirt. Clean it with a vacuum cleaner regularly.</p>								
Green	Quarterly	<p>Clean:</p> <p>The sealing strips around the freezer door must be cleaned regularly to prevent discolouration and prolong service life. Use clean water. After cleaning the sealing strips, check that it continues to provide a tight seal.</p>								

Maintenance chart

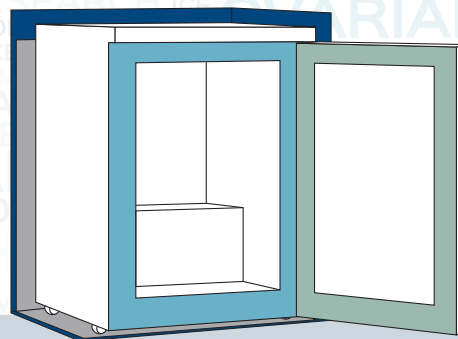

Maintenance chart

Dark blue

Weekly

Check placement:

Date:

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

Light blue

Weekly

Door openings

per 24h / Monthly

manual defrosting

< 5 / 1-2

< 10 / 2-4

< 15 / 4-6

Date:

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

Gray

Monthly

Clean:

Date:

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

Green

Quarterly

Clean:

Date:

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—

—/—/—