


# Maintenance for Chest Freezers

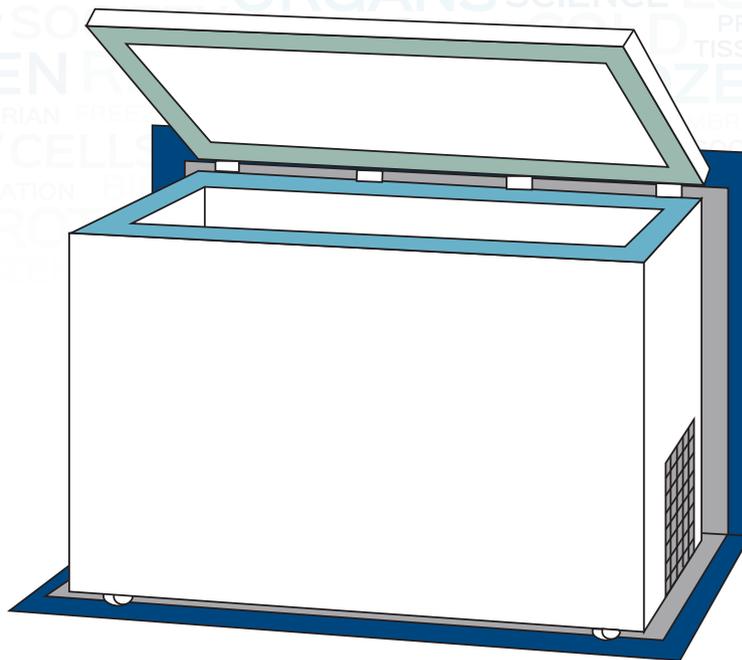

Maintenance										
Dark blue	Weekly	<p><b>Check placement:</b>                      Min.10 cm free space to the side of the freezer                      Min. 10 cm free space to the back of the freezer</p>								
Light blue	Weekly	<p><b>Defrosting:</b>                      Small amounts of loose ice can be removed using a plastic scraper.                      Defrosting has to be done regularly to prevent ice building                      The ice build depends on the use of the freezer.                      The more daily lid openings - the more manual defrosting.</p> <table border="0"> <tr> <td>Lid open/close per 24h</td> <td>Advised monthly manual defrosting</td> </tr> <tr> <td>&lt; 5</td> <td>1-2</td> </tr> <tr> <td>&lt; 10</td> <td>2-4</td> </tr> <tr> <td>&lt; 15</td> <td>4-6</td> </tr> </table>	Lid open/close per 24h	Advised monthly manual defrosting	< 5	1-2	< 10	2-4	< 15	4-6
Lid open/close per 24h	Advised monthly manual defrosting									
< 5	1-2									
< 10	2-4									
< 15	4-6									
Gray	Monthly	<p><b>Clean:</b>                      The ventilation grille, the condenser, the compressor, the compressor compartment and the floor must also be kept free of dust and dirt. Clean it with a vacuum cleaner regularly.</p>								
Green	Quarterly	<p><b>Clean:</b>                      The sealing strip around the freezer lid must be cleaned regularly to prevent discolouration and prolong service life. Use clean water. After cleaning the sealing strip, check that it continues to provide a tight seal.</p>								

